
CFH

No. 52175

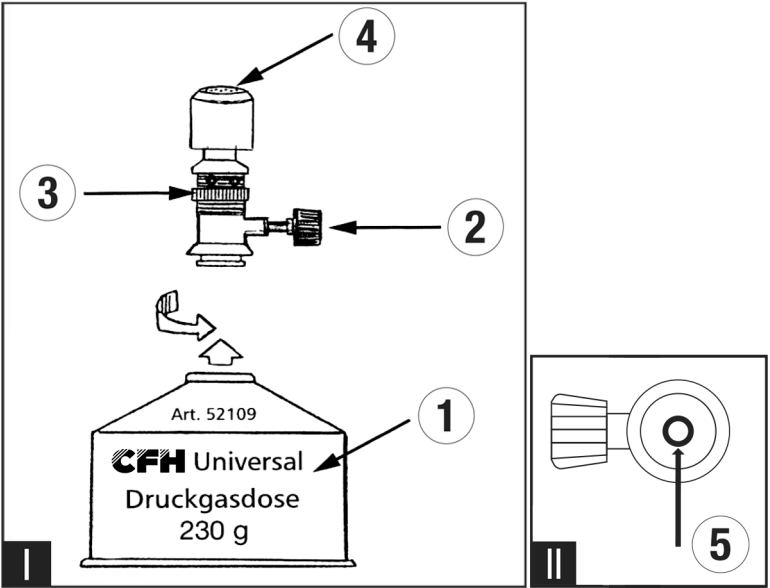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Bunsen-Laborbrenner BL 1700

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GB Operating manual

Bunsen Lab Burner BL 1700

Important! Read through this instruction manual carefully to become familiar with the device before you attach it to a gas canister. Keep the instructions in a safe place so that you can read them again.

Intended use

The Bunsen laboratory burner BL 1700 can be used universally and is ideal for the following work: Lead casting, goldsmith and laboratory work, test tube heating, optics, glass and plastic forming, physics and chemistry experiments, dental and pharmaceutical work.

This device may only be used with the 230 g (CFH part no. 52109) or 440 g (CFH part no. 52108) butane/propane gas cartridge. It could be dangerous to try to use other kinds of gas receptacles. This cartridge has an extraction valve with a connection thread that is standardised according to EN 417.

Operate the device in well-ventilated areas only.

Equipment

- (1) Pressurised gas container UG 230 (CFH part no. 52109)
- (2) Gas regulating valve for flame size adjustment
- (3) Air intake holes
- (4) Burner
- (5) Black rubber seal

Delivery

- 1 Bunsen laboratory burner
- 1 CFH pressurised gas container UG 230 (CFH part no. 52109)
- 1 User manual

Please check that the article is complete.

! Safety instructions

- Keep the Bunsen laboratory burner and the gas cartridge out of reach of children. The device must never be used by children (**danger of being injured**)!
- Operate in well ventilated spaces only. The use in closed rooms is forbidden.
- The gas canister is a disposable container and it cannot be refilled. Refilling is prohibited as it is very dangerous (**danger of being injured**).
- Always ensure that there is a seal fitted (between the device and the gas canister) before you attach the gas canister and that it is in a good condition (see Drawing II, Fig. 5 – the black rubber seal fitted on the device's gas extraction valve).
- The gas cartridge (No. 1) must always be firmly screwed onto the gas extraction valve of the device!
- Follow safety instructions stated on the cartridge.
- Ensure that the gas regulation valve (No. 2) is always tightly closed after you have finished working. Afterwards screw off the gas cartridge from the device.
- Do not make any changes to the gas cartridge or the device (**risk of injury**).
- Do not expose the gas canister to temperatures above 50°C (e.g. sunlight) (**danger of an explosion**).
- Only change the gas cartridge when outdoors and at a sufficient distance from other persons. Smoking is prohibited when changing canisters.
- The content of the gas cartridge is a mixture of propane/butane and is flammable. The device and the gas cartridge have to be protected from the reach of children and unauthorized persons.
- Never transport the gas cartridge in an airplane.
- Avoid using the device near flammable materials (paper, wood or inflammable substances). The device can reach very high temperatures during normal operation.
- CAUTION:** accessible parts can be very hot. Keep small children away from the device (**risk of burns**).
- Avoid touching the burner while it is hot.
- Do not put the ignited burner down (**danger of fire**). Do not leave the device unattended if it is on. Caution! Even after use, burners must never be placed near flammable materials – **danger of fire**.
- Never use the device in a barn, in a wooden gazebo or near a wooden fence (**risk of fire**).
- The device must be operated at a secure distance from flammable materials. A minimum distance of 50 cm from flammable substances must be observed.
- Do not use the device at all under conditions of strong wind, rain or high levels of humidity. If the flame is blown out by a gust of wind, immediately shut-off the gas regulating valve (No. 2), so that no more gas can flow out.
- Do not keep the device or the gas cartridge near heat sources (heaters, electrical appliances, ovens or strong sun radiation, etc).
- The device is only to be used in a well-ventilated area. National standards apply to:
 - supplying combustion air
 - to avoid the accumulation of dangerous amounts of unburnt gas.
- The device must never be used in the vicinity of heat sources, flames or sparks, dust accumulations or flammable substances. The device is only to be used at a safe distance from flammable materials – keep a safe distance of at least 0.5 m away from all objects as well as a distance of at least 1 m away from any objects that are above the device.

Symbol explanations

Open the gas regulating valve = Turn the gas regulating valve in the ((+ ON)) direction.

Close the gas regulating valve = Turn the gas regulating valve in the ((- OFF)) direction.

P = propane

Leak testing

If gas leaks out of your device (you can smell gas), take it out immediately to a place with a good air circulation with no inflammable sources, where you can look for the leak and fix it. Check the tightness of your device outdoors only. Never look for leaks with a flame. Use soapy water or a leak detection spray instead!

Important notes before putting the device into operation

Assembly and settings made by the manufacturer must never be altered. It is dangerous to change the structure of the device on your own authority, to remove parts or to use parts which are not approved by the manufacturer. Never modify the device in any way.

This device may only be operated with a gas cartridge with an extraction valve that complies with the European EN 417 standard (for example, the pressurised gas cartridge 230 g (CFH part no. 52109) or 440 g (CFH part no. 52108). It can be dangerous to try to connect it to other types of gas canisters. These cartridges can be obtained in building and specialized stores easily.

Do not use the device in a way where the gas cartridge is held horizontally or where the gas cartridge valve is pointing downwards. In this case, liquid gas would stream into the burner, which would lead to a reddish blazing uncontrolled flame. It is possible that the burner will then malfunction. The appliance has to be immediately positioned upright (see picture I) with the gas cartridge valve pointing up. At the same time the gas regulation valve (No. 2) must be closed. After 2-3 minutes the appliance will be fully functional again.

Do not shake the gas cartridge or the device with properly attached cartridge, before operation or during work. Shaking can cause the liquid gas to stream out. If this happens, close the gas regulating valve (No. 2) immediately. After 2-3 minutes the appliance is fully functional again, because the gas in the gas cartridge

has again been stabilized.

Change the gas container (cartridge) only when outdoors, far away from all possible combustible sources, as e.g. open fire, pilot lights, electric cooking appliances, and at a sufficient distance from other persons.

Please follow the safety instructions printed on the gas cartridge.

Measures to be taken before connecting the gas canister

Make sure that the terminals and connections are in place and damage-free. Before connecting the gas canister, make sure that the seal (no. 5 – drawing II) in the gas extraction valve of the device is present and in good condition (visual inspection). This gasket guarantees a gas tight connection between the device and the gas container (cartridge). Check the gas-carrying parts for damage. Do not use a device with a damaged or worn out gasket. Do not use a defective device that leaks or does not work properly. Before attaching the cartridge, close the gas regulation valve on the device. It is very important that always before you light the flame on the burner, you make sure that all parts are tightly connected. Never look for a leaking spot with a flame but use soapsuds (e.g. 10 drops of washing-up liquid with water) or a leak seeking spray.

Assembly of the pressurised gas container

The soldering torch is fully assembled except for the gas canister with its extraction valve. Make sure that you attach the gas cartridge (No. 1) to the device outside, away from possible combustible sources, as e.g. open fire, pilot flames, electrical cooking devices, and in a sufficient distance from other persons. Ensure that the gas regulating valve (No. 2) is tightly closed. Remove the protective cap from the gas canister before you attach it. A screwdriver is required, among other things. Now, holding the gas cartridge vertically, screw it TIGHTLY onto the gas extraction valve (see picture I). The rubber seal (no. 5) in the gas extraction valve must be checked before each new connection of the gas cartridge (drawing II – visual inspection). Check the tightness of the gas-carrying parts by brushing them with soapy water or spraying them with a leak-detection spray. Check the tightness of your device outdoors only. Never seek a leak with a flame, instead use soapy water or a leak detection spray (as described) for that!

Starting work

Close the gas regulation valve (No. 2). Remove the protective cover on the gas cartridge (No. 1). Then screw the gas cartridge (no. 1) TIGHTLY onto the gas extraction valve in the direction of the arrow. Then open the gas regulating valve (No. 2) slightly and ignite the gas at the outlet of the burner (No. 4) at the bottom with a lighter or a match. By turning the gas regulating valve (no. 2) in the (+) or (–) direction, you can determine the burner power setting. (IMPORTANT! Opening the gas regulating valve (No. 2) too wide will allow too much gas to flow into the burner (No. 4) and igniting the flame will then be difficult). After the flame has ignited, the gas regulating valve (No. 2) can be opened more to enable the device to reach full power. **Note: The cartridge is held by hand (this serves as a handle) during the intended use of the Bunsen laboratory burner. Always wear gloves if you are going to work with the soldering torch for a long time as the canister will cool down quickly in your hand while you are using it (up to 5 °C). When putting it down on a surface, you must ensure that you do not knock the device over, as the canister's base is not very wide.** The device must not be shaken before ignition and during preheating, as the gas may escape in the liquid phase. A preheating time of at least 3 minutes must be maintained until the working temperature is reached. Do not hold the burner above your head or tilt it down too much during preheating.

Especially important!

The burner must be preheated for at least 3 minutes until the burner head has reached the operating temperature. During the pre-heating period or due to strong shaking of the device the liquid gas can stream out.

Flame adjustment

On the one hand, you can regulate the gas flame intensity with the regulation valve (fig. 2). On the other hand, you also have the additional option of regulating the gas flame accordingly by turning the oxygen regulation ring (fig. 3). The hottest flame can be achieved with the air intake holes fully open. If you close these holes with the oxygen regulation ring (fig. 3), the flame intensity will be reduced.

Switching off

Turning the gas regulation valve (fig. 2) clockwise (rotation to the right) the gas supply stops and the flame extinguishes. Make sure in any case before taking off the gas cartridge that the flame at the burner has extinguished.

Replacing the gas container (gas canister with extraction valve)

Ensure the flame has been extinguished before you loosen the gas cartridge. Then unscrew the gas cartridge from the device and allow the device to cool completely before storing or transporting it. Check the seal (No. 5, Figure II - visual inspection) before connecting a new gas container. Change the gas container outdoors and away from other people. Before you attach the gas cartridge (No. 1) to the device - outside, far away from possible combustible sources, as e.g. open fire, pilot flames, electrical cooking devices, and in a sufficient distance from other people -, please ensure that the gas regulation valve (No. 2) is tightly closed. Now, holding the gas cartridge vertically, screw it tightly onto the gas extraction valve (No. 1), as shown above (picture I), so that no gas can stream out. The gasket must be checked before each new connection of the cartridge.

Cleaning the device

The device can be cleaned with a dry cloth when the burner is cool. Do not use any cleansing agents because the openings in the burner and the gas nozzle could become blocked.

Storage and transport

After end of work, before storage and transport, always unscrew the gas cartridge from the device. Important note: when you unscrew the gas canister, the canister's extraction valve will shut off automatically so that no more gas can escape

When the device is not used, store it protected from dust and humidity.

Ensure that the device has cooled down and the gas extraction valve is closed, before you store it securely and in a dry place. The same applies to transport. Make sure the device and the gas cartridge cannot move back and forth uncontrollably. Always transport the device and the gas cartridge unscrewed.

Maintenance + repair

Use the specified replacement parts only. Never operate a device with a damaged part. Any repairs as well as replacing the nozzle or replacing the gasket must only be carried out by an authorised specialist liquid gas dealer or by the manufacturer. Contact the manufacturer at the service address given below if repairs are needed.

Do not make any changes to the device. Assembly and settings made by the manufacturer must never be altered. It is dangerous to change the structure of the device on your own authority, to remove parts or to use parts which are not approved by the manufacturer.

Guarantee

Always keep the original receipt in a safe place. This document is needed as proof of purchase. Contact us via e-mail at: info@cfh-gmbh.de if you have a guarantee claim. We will contact you immediately.

Disposing of the gas canister with its extraction valve

A fully emptied pressurised gas canister can be taken to the collecting point in a dual recycling system for disposal. Partially emptied pressurised gas cans have to be disposed of at the collection centre for harmful substances.

Disposal of the device

The burner is made from recyclable materials. Contact your local recycling office or waste disposal company for disposing of or recycling the burner.

Technical specifications

Working temperature:	approx. 600 °C
Consumption:	approx. 145 g/h
Temperature threshold level:	approx. 2.0 kW
Type of gas:	Propane/butane mixture
Device category:	immediate pressure
Burner ø:	20 mm

The power rating during the normal operation can vary from the values set in the operating instructions (e.g. when using longer on large areas the performance declines. The reason for this is the icing up of the gas cartridge due to the withdrawal of gas). **Important information: The ambient air temperature will have a decisive impact on the performance of the device. In general, the following applies: the performance dissipates when it is cold.**

We reserve the right to make technical and visual changes.

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